# Explorer<sup>®</sup> Semi-Micro Balances

## User Guide

EN

ES

FR

DE

IT

PT

SV

NL

DA

PL

CZ

HU

## 1. SAFETY INFORMATION

**WARNING** For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.

**CAUTION** For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.

**Note** For useful information about the product

General Hazard

Electric Shock Hazard

### Safety Precautions

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

#### Safety precautions for the Ionizer:

**WARNING: ELECTRICAL SHOCK HAZARD**  
AVOID TOUCHING THE ELECTRODE PINS WHILE THE IONIZER IS TURNED ON.

#### Intended Use

Use the instrument exclusively for weighing determination as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

## 2. INSTALLATION

### 2.1 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes. Allow sufficient space around the instrument.

## 2.2 Installing the Wind Ring, Weighing Platform

## 2.3 Installing the Draft Shield Glass and Doors

### Side Door Installation

1. Unlock
2. Align
3. Clear
4. Lock

### Front Panel Installation

1. Release terminal
2. Align
3. Push to rear
4. Reconnect terminal


### Top Door Installation

1. Align
2. Press Down

## 2.4 Installing the Glass Shelf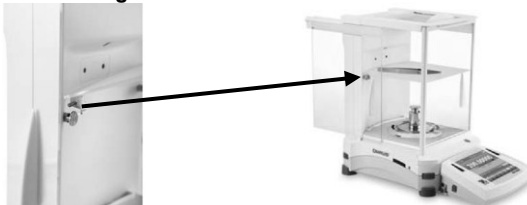

## 2.5 Connecting Power

### Models with AC Adapter

Connect the AC adapter power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

**Attention:** Only use an AC adapter specified by OHAUS.


**Note:** Allow equipment to warm up for at least 2 hours for optimal weighing performance.

## 2.6 Leveling the Equipment

To level the balance, adjust the **Leveling Thumbwheels** at each corner until the bubble is centered in the circle.

Be sure the equipment is level each time its location is changed.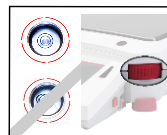

## 2.7 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

# 3. OPERATION

## 3.1 Overview of Display, Home Screen

This equipment utilizes a touch-sensitive display. *Touch* areas and Buttons to control the equipment's functions.

### CONTROLS

#### Instructional Messages

Stability (\*), Net (NET), Gross (G) and/or center of zero (>0<) indicators

Result Fields: Information varies by application

Touch icons to perform specific functions or access other functions

Touch to access application setup menu

Touch to change applications

Touch to access available libraries

Dynamic Weighing Library

Place sample on the pan. Press Begin Dynamic to start averaging.

\* 30.07477 g

Capacity Guide 0 220

Averaging Time: 10 s

Edit Settings Tare Begin Dynamic

Print Zero Sensors Ionizer AutoCal™ Menu More...

Touch for information menu

Touch unit indicator to change weighing unit

Capacity or Status Guides: Vary by application

Application Buttons: Functions vary by application

## 3.2 Principal Functions and Main Menu

**Weighing:** Press **Tare** to set the display to zero. Place an item on the pan. Display indicates gross weight.

**Taring:** With no load on the pan, press **Tare** to set the display to zero. Place an empty container on the pan and press **Tare**. Add material to the container and its net weight is displayed. Remove container and container's weight appears as a negative number. Press **Tare** to clear.

## MENU & SCREEN NAVIGATION

Touch **Menu** to open the menu list.  
 Touch and drag the **Scroll bar** to view additional items.

**Calibration:**  
 Touch to view calibration options.

**User Settings:**  
 Touch to view user preferences.

**Balance Setup:**  
 Touch to view balance settings.

**Application Modes:**  
 Touch to view application modes.

**Weighing Units:**  
 Touch to view weighing units.

**GLP and GMP Data:**  
 Insert user data for traceability.

**Communication:**  
 Touch to view COM Device Settings and Print Settings.

**Library:**  
 Touch to delete all Library records.

**I/O Settings:**  
 Touch to view remote I/O Settings.

**Diagnostics:**  
 Touch to view Diagnostics items.

**Factory Reset:**  
 Touch to do a Factory reset.

**Lockout:**  
 Touch to access the Lockout menu list.

**Information:**  
 Touch to view list of topics.

### 3.3 Using the Balance

#### Weighing Application

1. If required, place an empty container on the pan and press **Tare**.
2. Place the object to be weighed on pan (or in container) to display the weight. When stable, the \* appears.
3. The resulting value is displayed in the active unit.

## 4. MAINTENANCE

### 4.1 Cleaning

**WARNING:** Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

**Attention:** Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

### 4.2 Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website [www.ohaus.com](http://www.ohaus.com) to locate the Ohaus office nearest you.

## 5. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only
- Altitude: Up to 2000 m
- Operating Temperature range: 5°C to 40°C
- Specified Temperature range: 15°C to 25°C
- Humidity: maximum relative humidity 80 % for temperatures up to 30°C decreasing linearly to 50% relative humidity at 40°C
- Mains supply voltage fluctuations: up to ±10% of the nominal voltage
- Installation category II
- Pollution degree: 2

**SPECIFICATIONS :**

MODEL	EX125D	EX125	EX225D	EX225D/AD	EX225/AD
Capacity	52g/120g	120g	120g/220g	120g/220g	220g
Readability d	0.01mg/0.1mg	0.01mg	0.01mg/0.1mg	0.01mg/0.1mg	0.01mg
Repeatability (std. dev.) (20g)	0.015mg	0.015mg	0.015mg	0.015mg	0.015mg
Repeatability (std. dev.) (100g)	0.1mg/0.1mg	0.02mg	0.02mg/0.1mg	0.02mg/0.1mg	0.03mg
Linearity (g)	±0.1mg	±0.1mg	±0.1mg	±0.1mg	±0.1mg
Power supply	AC Adapter Input: 100-240 VAC 0.6A 50-60 Hz AC Adapter Output: 12 VDC 1.5A				

MODEL	EX125DM	EX125M	EX225DM	EX225DM/AD	EX225M/AD
Capacity	52g/120g	120g	120g/220g	120g/220g	220g
Readability d	0.01mg/0.1mg	0.01mg	0.01mg/0.1mg	0.01mg/0.1mg	0.01mg
Verification scale interval, e	1mg	1mg	1mg	1mg	1mg
Repeatability (std. dev.) (20g)	0.015mg	0.015mg	0.015mg	0.015mg	0.015mg
Repeatability (std. dev.) (100g)	0.1mg/0.1mg	0.02mg	0.02mg/0.1mg	0.02mg/0.1mg	0.03mg
Linearity (g)	±0.1mg	±0.1mg	±0.1mg	±0.1mg	±0.1mg
Power supply	AC Adapter Input: 100-240 VAC 0.6A 50-60 Hz AC Adapter Output: 12 VDC 1.5A				

**6. COMPLIANCE**

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at <a href="http://www.ohaus.com/ce">www.ohaus.com/ce</a> .
	This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to <a href="http://www.ohaus.com/weee">www.ohaus.com/weee</a> .

**Verified weighing instruments**

When the instrument is used in trade or a legally controlled application it must be set up, verified and sealed in accordance with local weights and measures regulations. It is the responsibility of the purchaser to ensure that all pertinent legal requirements are met.

Weighing Instruments verified at the place of manufacture bear the following supplementary metrology marking on the descriptive plate.

Weighing Instruments to be verified in two stages have no supplementary metrology marking on the descriptive plate. The second stage of conformity assessment must be carried out by the applicable weights and measures authorities. If national regulations limit the validity period of the verification, the user of the weighing instrument must strictly observe the re-verification period and inform the weights and measures authorities. As verification requirements vary by jurisdiction, the purchaser should contact their local weights and measures office if they are not familiar with the requirements.

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P/N 30283440 A © 2015 Ohaus Corporation, all rights reserved.
Ohaus Corporation